TE70X Semi-dry Blotters

TE70X Semi-dry transfer unit
TE77X Semi-dry transfer unit
TE70XP Semi-dry transfer unit
TE77XP Semi-dry transfer unit

Contents

| | |
|---|-----|
| Important Information | ii |
| Waste Electrical and Electronic Equipment (WEEE) | vii |
| 1. TE70X Semi-dry Blotters: Description | 1 |
| Unpacking | 2 |
| Specifications | 3 |
| 2. Operating instructions | 4 |
| Electrotransfer | 9 |
| 3. Care and maintenance | 13 |
| 4. Troubleshooting | 14 |
| 5. Electrotransfer notes..... | 17 |
| 6. Bibliography..... | 19 |
| 7. Customer service information..... | 20 |
| 8. Ordering information..... | 21 |

Important Information – English

- If this equipment is used in a manner not specified by Hoefer, Inc. the protection provided by the equipment may be impaired.
- This instrument is designed for indoor laboratory use only.
- Only accessories and parts approved or supplied by Hoefer, Inc. may be used for operating, maintaining, and servicing this product.
- **Warning!** Because this instrument can develop sufficient voltage and current to produce a lethal shock, care must be exercised in its operation.
- This instrument is designed in accordance with the EN61010-1:2001 electrical safety standard. Nevertheless, it should be used only by properly trained operators. Read this entire manual before using the instrument and use only according to the instructions.
- The instrument must always be used with the earth lead of the power cord correctly grounded to earth at the mains outlet.
- Use only undamaged electrical wire and equipment specific for the voltages you will use. All equipment connected to high voltage should be in accordance with EN61010-1:2001.
- Keep the instrument as dry and clean as possible. Wipe regularly with a soft, damp cloth. Let the instrument dry completely before use.
- Do not operate the instrument in extreme humidity (above 80%). Avoid condensation by letting the unit equilibrate to ambient temperature when taking the instrument from a colder to a warmer environment.
- To permit sufficient cooling, ensure that the vents of the instrument are not covered.

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Waste Electrical and Electronic Equipment (WEEE)

English

This symbol indicates that the waste of electrical and electronic equipment must not be disposed as unsorted municipal waste and must be collected separately. Please contact an authorized representative of the manufacturer for information concerning the decommissioning of your equipment.

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
TE70X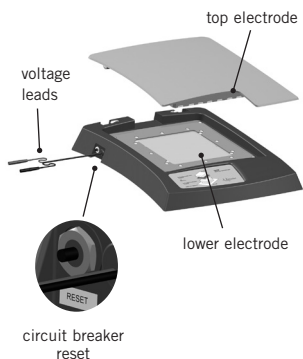

TE77XP

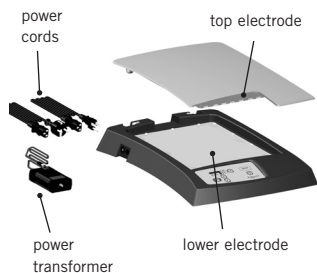
Fig 1. Semi-dry transfer unit main components.

The TE70X, TE77X, TE70XP, and the TE77XP semi-dry blotters rapidly transfer proteins from polyacrylamide gels onto a membrane by means of a low current and low voltage electrotransfer with minimal Joule heating. Most transfers are complete in one hour or less.

The smaller TE70X unit transfer surface is 14 × 16 cm, suitable for transferring standard gels, including those from the SE600 Chroma and the SE400. The larger TE77X transfer surface is 21 × 26 cm, suitable for transferring large format gels.

The TE70X and TE77X have safety circuits built into the instrument that limit the voltage to 30 V and the current to 0.5 A. The circuits protect the user from unnecessary electrical hazards.

The TE70XP and the TE77XP have a built-in power supply for transferring gels. This eliminates the need for an external power supply. These instruments can deliver up to 30 V and 0.5 A. The instruments also monitor the transfer stack resistance, and can stop a transfer if large changes indicate that the buffer system is depleting.

Multiple gels can be transferred at the same time by placing several small gels of the same thickness side by side, or by stacking two gels vertically in a carefully constructed multi-layered stack.

The electrodes are made of the best possible materials to ensure the units last a long time.

Unpacking

Unwrap all packages carefully and compare contents with the packing list, making sure all items arrived. If any part is missing, contact Hoefler, Inc. Inspect all components for damage that may have occurred while the unit was in transit. If any part appears damaged, contact the carrier immediately. Be sure to keep all packing material for damage claims or to use should it become necessary to return the unit.

Specifications

This declaration of conformity is only valid for the instrument when it is:

- used in laboratory locations.
- used as delivered from Hofer, Inc. except for alterations described in the user manual.
- connected to other CE labeled instruments or products recommended or approved by Hofer, Inc.

| | TE70X | TE77X | TE70XP | TE77XP |
|---------------------|----------|----------|--------------|--------------|
| Electrode size (cm) | 14 × 16 | 21 × 26 | 14 × 16 | 21 × 26 |
| <i>Inputs:</i> | | | | |
| Voltage (V) | 30 (DC)* | 30 (DC)* | 100–240 (AC) | 100–240 (AC) |
| Current (A) | 0.5* | 0.5* | 0.7 | 0.7 |
| <i>Output:</i> | | | | |
| Voltage (VDC) | 30 | 30 | 30 | 30 |
| Current (A) | 0.5 | 0.5 | 0.5 | 0.5 |
| Frequency (Hz) | | | 47–63 | 47–63 |

*Maximum setting for the customer's power supply.

Minimum transfer stack thickness:

| | |
|---------------|------------------|
| TE70X, TE70XP | 3.2 mm / 12.0 mm |
| TE77X, TE77XP | 2.8 mm / 12.0 mm |

| | |
|-------------------------------------|--|
| Environmental operating conditions: | Indoor use: 4–40 °C Humidity up to: 80% Altitude up to: 2000 m Installation category: II Pollution degree: 2 |
|-------------------------------------|--|

| | |
|-------------------------|--------------------------------------|
| Dimensions (w × d × h): | 38 × 46 × 9 cm (15 × 18 × 3.5 in) |
|-------------------------|--------------------------------------|

| | |
|---------|--------------------------------|
| Weight: | Shipping 6.8 kg Unit 3.7 kg |
|---------|--------------------------------|

| | |
|----------------|--|
| Certification: | EN61010-1, EN 61326, CE UL61010-1-2004 CSA 22.2 61010-1-04 |
|----------------|--|

2. Operating instructions

To transfer proteins, prepare the unit, assemble the stack, and connect to a power supply if necessary. Then run the transfer for the required amount of time. Each step is described below.

①

Prepare the unit by rinsing the electrodes with distilled water.

②

Prepare the gel

Cut away the wells and/or stacking gel section. Equilibrate the gel in transfer buffer if required.

③

Prepare the transfer stack

Cut the blot paper and transfer membrane to the same size as the gel. Stack the layers carefully so the edges align. If for some reason the membrane needs to be larger than the gel, use a mylar mask (see optional step below) to ensure the current does not bypass the gel.

Optional: Cut a mask to the proper size

Measure the gel. Cut an opening centered in the solid mask to a size roughly 2 mm smaller than the gel. Take care with sharp blades when cutting the masks. Place the mask in the base of the unit, centering the opening.

Note: Take care to place the gel correctly on the first try because proteins begin to transfer immediately; once transfer has begun, moving the gel will distort results or cause “shadow” bands on the blot.

4

Prepare the blotting paper

For each gel, cut at least 6 pieces of blotting paper the *same size as the gel* or *slightly smaller*.

Gauge the thickness or number of blotting paper layers according to the amount of buffer required; up to 300 ml of buffer may be required for larger gels or for transfers of 60 minutes in order to prevent the stack from drying out or the buffer from becoming depleted.

Saturate at least 3 pieces of the blotting paper with transfer buffer. One by one, center each sheet on the lower electrode and remove all trapped air by rolling a clean pipet or roller from the center toward the edges.

5

Prepare the membrane

For each gel, cut 1 membrane the same size as the gel or slightly smaller. (A larger membrane may contact an electrode panel, creating a pathway by which current can bypass the gel.)

Pre-wet nitrocellulose or nylon membranes with distilled water. Pre-wet PVDF or other hydrophobic membranes with methanol. Then soak all membrane types in transfer buffer for 2–5 minutes.

Note: Each 21 × 26 cm blot paper will absorb approximately 50 ml of transfer buffer. Each 14 × 16 cm blot paper will absorb approximately 20 ml of buffer.

Note: If not using Hoefer blotting paper, additional pieces may be required to achieve desired results.

Note: Always wear gloves when handling membranes to avoid leaving fingerprints.

6


Important! Stack each layer with care, with edges parallel. As each layer is added, remove all air pockets by rolling a clean pipet from the center to the edges. Add a few drops of buffer to trouble areas to help remove air pockets.

Note: If not using Hoefer blotting paper, additional pieces may be required to achieve desired results.

Complete the stack

- Place the pre-wet membrane onto the stack of blotting paper.
- Place the gel on the membrane.
Note: Proteins bind to the membrane as soon as contact occurs, so it is important to place the gel correctly on the first try.
- Cover the gel with three layers of buffer saturated blotting paper.

Fig 2. Transfer stack for a single gel.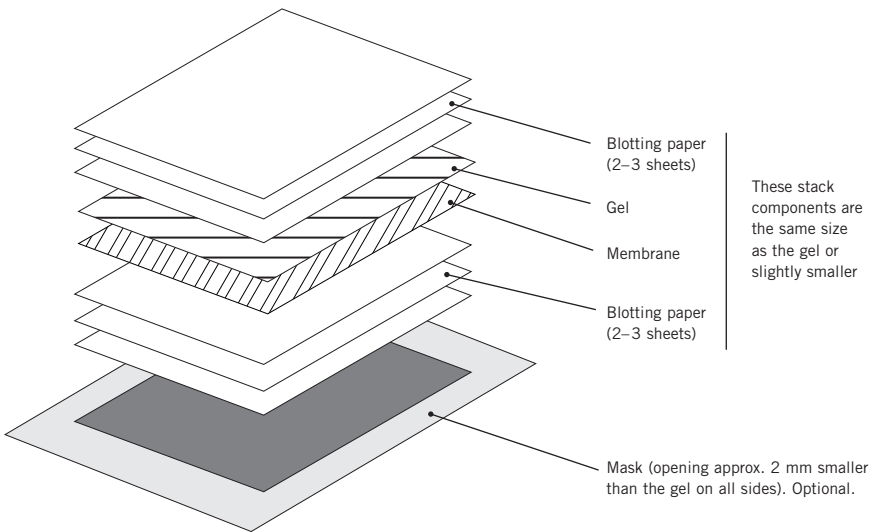

Note: When transferring multiple gels, transfer efficiency depends on such factors as gel thickness, gel position in the stack, transfer buffer, membrane type, and, most importantly, the characteristics of the protein. The gel closest to the anode generally transfers the most completely. It is preferable to lay gels side-by-side rather than stacking them.

Multiple gels: Either lay gels of the same thickness side-by-side (Fig 3), or stack 2 sandwiches layered as shown (Fig 4).

For best results, the transfer stack should be centered in the electrode panels.

If two gels are stacked, separate them with porous cellophane — **not plastic wrap!** (Cellophane permits electric current to pass but stops proteins.) Cut the cellophane slightly smaller than the gel and wet with transfer buffer.

Several sheets of buffer-soaked blotting paper on *each* gel provides electrical continuity.

7

Place the Cover on top of the transfer stack. Do not remove the cover until after the transfer is complete in order to prevent stack components from moving.

Fig 3. Transfer stacks for gels placed side-by-side.

Note: There must be no electrical contact between the two stacks.


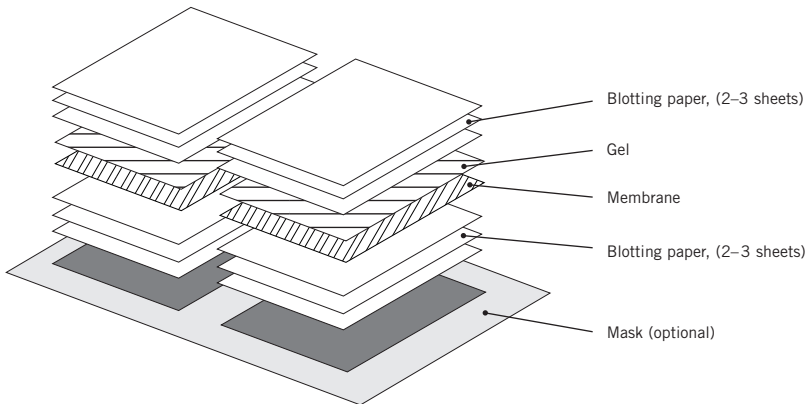
Fig 4. Transfer stack for stacked gels.

Separate the stack for each gel with a sheet of cellophane.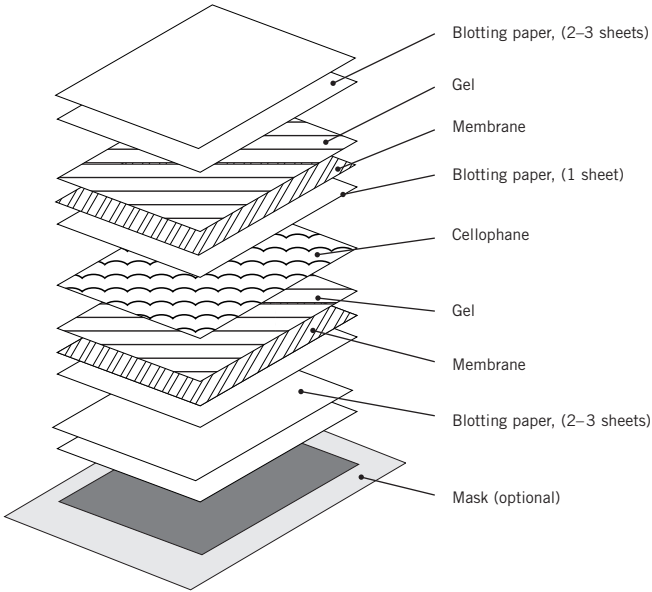

Electrotransfer

TE70X and TE77X

1

The power supply should be switched off and both the current and voltage controls set at zero

Then plug the color-coded leads from the base of the transfer unit into the power supply jacks, matching red to red and black to black. Do not reverse polarity. The lid contains the black, or negative electrode. The base contains the red, or positive electrode.

2

Set the power supply voltage and current

Set the voltage limit to 30 V maximum. The input voltage must not exceed 30 V or the protection will be activated. The maximum current should not exceed 0.8 mA/cm² of the gel surface or 500 mA maximum. If transferring several layers of gels, a longer time may be required.

3

If available, set the power supply timer


Most transfers are complete within one hour, but larger proteins, proteins from native gels, and thicker gels may require an additional 1 hour of transfer time. The optimum transfer time for each protein and gel system must be determined empirically.

Note: Transfers exceeding one hour may require additional sheets of buffer saturated blot paper in the transfer stack.

Note: Generally, smaller fragments transfer more quickly than larger ones.

Fig 5. Recommended current settings for different gel sizes.

Use the graph to quickly find the current setting for your gel size, or calculate the gel area (cm²) and multiply it by 0.8 mA.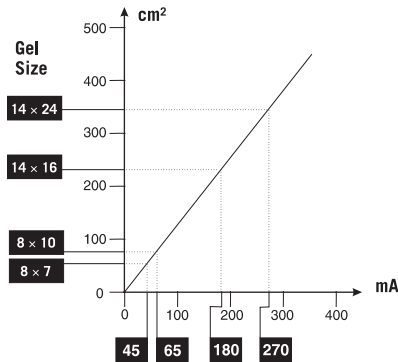

TE70XP and TE77XP

1

Note: While setting the time, there is an extra setting labelled "ON" above the 4:00 upper limit. This will disable the timer, allowing the transfer to run continuously until manually turned off by the user.

Turn the instrument on using the POWER button on the front keypad.

The instrument works by setting the current and time, and then starting the transfer. During the transfer, the voltage can be displayed. The transfer cannot be set to run at constant voltage.

The DISPLAY MODE BUTTON toggles between the current (mA), time (hour:minutes) and the voltage (V).


The UP and DOWN arrows to change the value for the current and time.

2

Press the START/STOP button to start the transfer.

| Variable | Units | Range | Increment |
|----------|---------------|--------------------|-----------|
| Current | milliamps | 1–500 | 1 mA |
| Time | hours:minutes | 5 min – 4:00 hours | 5 min |
| Voltage | volts | 0–30 (read only) | 1 volt |


Fig 6. TE70XP keypad.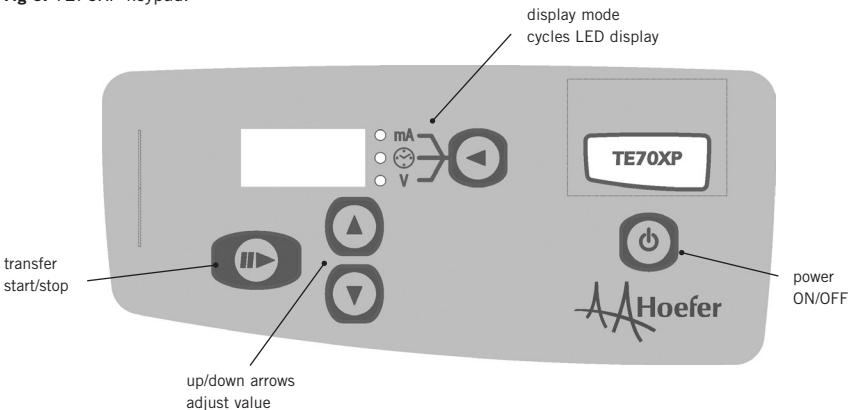

During the transfer

TE70XP and the TE77XP

A red LED will appear in the lower right corner of the Display when voltage is applied to the transfer stack.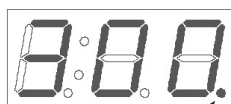

Red LED

Both the Time and the Current can be changed as the transfer progresses. Select mA or time using the DISPLAY MODE button, and the UP and DOWN to change the value. While the value is being changed, the LED will blink. After about 10 seconds the LED will stop flashing and show the real time run parameters.


Buffer Depletion

One of the more common failure modes for a Western transfer is buffer depletion. Buffer depletion leads to changes in pH and overheating, both of which are detrimental to the transfer.

The TE70XP and the TE77XP instruments monitor the transfer stack resistance. Large changes in this resistance indicate the depletion of the buffer system. The instrument can stop a transfer before these changes lead to further problems, like burning of the transfer stack.

If this condition happens, the error message “dry” will be on the display. The instrument has stopped the transfer. If desired, the cover can be removed, and buffer can be added to the transfer stack. Replace cover, and press any key (except Power ON/OFF) to continue.

For future transfers, add more layers of buffer soaked blot paper.

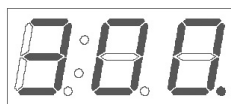Dry error message.

After the transfer is complete

TE70X and TE77X

Turn off the power supply and disconnect the leads from the power supply.

TE70XP and TE77XP

After a transfer is complete, the unit will beep for 5 seconds. The LED display will blink and cycle between the values for the current, time and voltage at the time the transfer was stopped. These can be recorded into a notebook, if desired. Pushing any button will clear the display.

1

Remove the cover slowly because the stack may adhere to it.

2

Remove and dispose of the upper blot papers.

3

Remove the gel(s).

Optional: Stain gel to check for residual protein left in the gel.

4

If desired, label the gel contact side of the membrane with a soft pencil.

5

Remove the membrane(s) from the stack with blunt forceps. Process the membrane according to your protocol or allow the membrane to air dry prior to storage.

6

Remove the remaining blot papers and dispose.

7

Rinse the unit according to the Care and maintenance instructions (next page).

Note: Staining the gel(s) for residual protein gives an indication of the completeness of transfer.

Note: Rewet dried membranes in the appropriate wetting buffer prior to processing.

3. Care and maintenance

- Do not autoclave or wash the unit in a dishwasher.
- Do not immerse the unit in water!

Rinse the cathode in the cover and the anode in the base with distilled water. Let the unit air dry completely. If using radioactive reagents, decontaminate the unit with a cleaning agent such as Contrad® 70. Never use abrasive cleansers.

4. Troubleshooting

problem

solution

Incomplete transfer

Blank or faint areas on the membrane

Remove trapped air pockets between the gel and membrane during stack assembly.

Use buffer with a lower ionic strength.

Molecules do not migrate out of gel

Check all electrical connections. Confirm that current is flowing through the transfer stack.

Check that the buffer pH is close to the intended pH. Most buffers should not be titrated. Make fresh buffer.

Use 3.5 mM SDS (0.1%) in the transfer buffer.

Add several more sheets of buffer-saturated blotting paper to each side of the gel sandwich so that more buffer is present during the transfer.

Increase the transfer period. Large fragments may require an additional hour.

Do not use staining or fixing agents on the gel before transfer.

Use a thinner gel.

Reduce the gel acrylamide concentration.

If using a non-nitrocellulose membrane, avoid including methanol in the transfer buffer or reduce the amount to the minimum possible.

Use reagent-grade chemicals.

Increase the net charge on the protein by using a transfer buffer with a different pH. Lower pH (<6–7) increases the positive charge on proteins; higher pH (>6–7) increases the negative charge on proteins.

problem**solution**

Open circuit or no output current for TE70X or TE77X

Push circuit protection reset button located next to the input leads. To prevent recurrence limit the input voltage to 30 V and the current to 500 mA.

Early models: Blown fuse in protection circuit. Replace 1.6 A micro fuse. (See Accessories in Ordering information.)

Smearred or diffuse band patterns

If equilibrating before the transfer, shorten or eliminate the equilibration time and/or equilibrate under coldroom conditions.

If the transfer buffer contains $\geq 10\%$ methanol, equilibrate the gel in transfer buffer for 30 minutes to allow it to shrink before assembling the stack.

Note: Large proteins may not migrate as readily once the pore size is slightly reduced.

Take care that the gel does not shift once it contacts the membrane.

Buffer depletion can change the pH in the transfer stack, and have a negative effect on the transfer. On subsequent transfers, either shorten the transfer time, reduce the current, or increase the number of buffer-soaked blotting papers in the stack.

Check that the preferred binding surface of the membrane (if any) contacts the gel.

Uneven band transfer

The blotting paper and membrane must be the same size as the gel or 1–2 mm smaller. Larger sizes will provide an electrical path for current to bypass the gel solution.

Different proteins will transfer at different rates depending on size and net charge.

problem**solution**

Inefficient binding to membrane*Chemical parameters*

Prepare transfer buffer without SDS. (SDS can improve transfer efficiency but can interfere with protein binding to a PVDF membrane.)

Add 10–20% methanol to the transfer buffer to enhance binding to nitrocellulose or PVDF.

Membrane parameters

Use a membrane with a smaller pore size (0.20 μm) if proteins pass through the membrane.

Place a membrane both over and under the gel to capture any proteins migrating in the opposite direction.

Check if too much sample is available for the binding surface area by placing two membranes instead of one. If “blow through” occurs, reduce the sample load.

Wear gloves when handling membranes.

Store membranes at ambient temperature and out of direct sunlight.

Check shelf life of nitrocellulose membrane, replace if necessary.

TE70XP, TE77XP*Current reading lower than setting*

Maximum voltage (30 V) has been reached.
Buffer may be depleted.

No output current

Transfer stack is not making contact with the upper electrode. Add buffer soaked blot papers.

“dry” error message

Instrument has detected large changes in transfer stack resistance.

Add more buffer soaked blot papers in future transfers.

No output voltage

Transfer stack is being short circuited.

Note: For more troubleshooting ideas, refer to Bjerrum, O.J. *et al.* (1988).

5. Electrotransfer notes

- Run the transfer as soon as possible after electrophoresis to minimize protein diffusion within the gel.
- Stacked gels must all be the same size.
- Limit transfers to two hours or less.
- The recommended methanol concentration for different membrane types are:

| membrane type | methanol % |
|----------------|------------|
| Charged nylon | 0 |
| Nitrocellulose | 10–20 |
| PVDF | 10–20 |

- Use a buffer with low ionic strength such as one of the two listed below to prevent overheating. Use the CAPS buffer when Tris cannot be used (*e.g.*, peptide sequencing). CAPS can improve transfer because of its effect on the charge of the protein (see Matsudaira, 1987).

Towbin buffer

(25 mM Tris, 192 mM glycine, 20% v/v methanol, pH 8.3, 1 liter)

| | | |
|--------------------|---------------|--------|
| Tris (FW 121.1) | 25 mM | 3.0 g |
| Glycine (FW 75.07) | 192 mM | 14.4 g |
| SDS* (FW 288.4) | 0.1% (3.5 mM) | 1.0 g |

Dissolve in 600 ml distilled water.

Add methanol as required[†].

Bring to 1 liter with distilled water. Do not adjust the pH, which should be between 8.2–8.4.

Optional: Chill before use.

**Optional:* Adding SDS can improve transfer efficiency.

[†]Depending on the membrane type selected (see table above), adding methanol can improve transfer results.

Note: Buffers containing methanol may deteriorate if stored for long periods — add methanol just prior to transfer.

CAPS buffer, 1X

(10 mM CAPS, pH 11.0, 1 liter)

| | | |
|--|-------|-------|
| CAPS (FW 221.3) | 10 mM | 2.2 g |
| [3-(cyclohexylamino)-1-propanesulfonic acid] | | |

Dissolve in 600 ml distilled water, adjust to pH 11.0 with conc. NaOH.

Adjust volume to 1.0 liter.

- For a 3-buffer system, refer to Kyhse-Anderson, J. (1984).
- Transfer efficiency varies depending on the gel concentration, which can be optimized. For more information, refer to Smejkal and Gallagher (1994).

6. Bibliography

- Bjerrum, O.J., Larsen, K., and Heegaard, N., *CRC Handbook of Immunoblotting of Proteins 1*, Section 7. CRC Press (1988).
- Gallagher, S., Winston, S.E., Fuller, S.A. and Hurrell, J.G.R., Immunoblotting and Immunodetection. In *Current Protocols in Molecular Biology*. 10.8.1–10.8.17. Greene Publishing and Wiley-Interscience, NY (1993).
- Hancock, K. and Tsang, V., India ink staining of proteins on nitrocellulose paper. *Anal. Biochem.* **133**, 157–162 (1983).
- Kyhse-Anderson, J., Electroblotting of multiple gels: A simple apparatus without buffer tank for rapid transfer of proteins from polyacrylamide to nitrocellulose. *J. Biochem. and Biophys. Meth.* **10**, 203–209 (1984).
- Matsudaira, P., Sequence from Picomole Quantities of Proteins Electroblotted onto Polyvinylidene Difluoride Membranes. *J. Biol. Chem.* **262**, 10035 (1987).
- Sasse, J. and Gallagher, S., Detection of Proteins on Blot Transfer Membranes. In *Current Protocols in Molecular Biology*. 10.7.1–10.7.3. Greene Publishing and Wiley-Interscience, NY (1991).
- Smejkal, G., and Gallagher, S., Determination of Semidry Protein Transfer Efficiency with Transverse Gradient Gel Electrophoresis. *Biotechniques*. **16**, 196–202 (1994).
- Tovey, E. and Baldo, B., Comparison of semi-dry and conventional tank-buffer electrotransfer of proteins from polyacrylamide gels to nitrocellulose membranes. *Electrophoresis* **8**, 384–387 (1987).

7. Customer service information

Technical service and repair

Hoefler, Inc. offers complete technical support for all of our products. If you have any questions about how to use this product, or would like to arrange to repair it, please call or fax your local Hoefler, Inc. representative.

Check the Hoefler, Inc. website at www.hoeflerinc.com for the distributor in your area. Or contact us directly at:

Hoefler, Inc.

84 October Hill Road
Holliston, MA 01746

Toll Free: 1-800-227-4750

Phone: 1-508-893-8999

Fax: 1-508-893-0176

support@hoeflerinc.com

www.hoeflerinc.com

Important! Request a copy of the Hoefler, Inc. "Health and Safety Declaration" form before returning the item. No items can be accepted for servicing or return unless this form is properly completed.

Note: A Return Authorization (RA) number must be obtained from Hoefler, Inc. before returning any item to Hoefler, Inc.

8. Ordering information

| product | qty. | code no. |
|--|------|----------|
| TE70X Semi-dry Transfer Unit, 14 × 16 cm Includes 25 sheets of blotting paper, 50 sheets of cellophane and 2 solid masks. | 1 | TE70X |
| TE77X Semi-dry Transfer Unit, 21 × 26 cm Includes 25 sheets of blotting paper, 50 sheets of cellophane and 2 solid masks. | 1 | TE77X |
| TE70XP Semi-dry Transfer Unit, 14 × 16 cm Includes 25 sheets of blotting paper, 50 sheets of cellophane and 2 solid masks. | 1 | TE70XP |
| TE77XP Semi-dry Transfer Unit, 21 × 26 cm Includes 25 sheets of blotting paper, 50 sheets of cellophane and 2 solid masks. | 1 | TE77XP |

Accessories

TE70X and TE70XP

| | | |
|------------------------------------|----|-----------|
| Solid masks, 16.5 × 18.5 cm | 4 | TE74 |
| Porous cellophane, 20 × 35.5 cm | 50 | TE73 |
| Blotting paper, precut, 14 × 16 cm | 25 | TE76-1416 |

TE77X and TE77XP

| | | |
|------------------------------------|----|--------|
| Solid masks, 23 × 27.5 cm | 4 | TE78 |
| Porous cellophane, 35 × 44 cm | 50 | SE1142 |
| Blotting paper, precut, 21 × 26 cm | 25 | TE76 |

| | | |
|---------------------------|-------|---------------|
| Power Cord Kit | 1 set | PS36-24 |
| Fuse, Micro, 1.6 A, 250 V | 5 | PSF1.6A-MICRO |

Related products

| product | code no. |
|---|---------------|
| SE600 Chroma | SE600X-15-1.5 |
| MiniVE Vertical Electrophoresis system | SE300-10A-1.0 |
| PBS300B Power Supply, 300 V, 500 mA, 90 W | PBS300B |

Transfer Membranes

Pure Nitrocellulose, sheets and roll

0.45 µm pore size

| | |
|-------------------------|--------------|
| 8 × 9.5 cm, 10 sheets | GM-NC45-89 |
| 16 × 16 cm, 10 sheets | GM-NC45-1616 |
| 20 × 23.5 cm, 10 sheets | GM-NC45-2320 |
| 30 cm × 3 m, 1 roll | GM-NC45 |

0.2 µm pore size

| | |
|---------------------|---------|
| 30 cm × 3 m, 1 roll | GM-NC22 |
|---------------------|---------|

PVDF membrane

0.45 µm pore size

| | |
|-----------------------|--------------|
| 10 × 10 cm, 10 sheets | GM-PV45-1010 |
| 15 × 15 cm, 5 sheets | GM-PV45-1515 |
| 30 cm × 3 m, 1 roll | GM-PV45 |

Hoefer Electrophoresis Reagents

| | |
|--|-----------|
| Tris, 1 kg | GR132-1 |
| Glycine, 1 kg | GR125-1 |
| SDS, 500 g | GR126-500 |
| Tween® 20, 500 ml | GR128-500 |
| DTT, 5 g | GR122-5 |
| Glycerol, 1 L | GR124-1 |
| Bromophenol Blue (BPB), 10 g | GR120-10 |
| Protein Determination Reagent, 500 standard assays | GR133-500 |
| Coomassie® Brilliant Blue G-250, 25 g | GR134-25 |
| Coomassie Brilliant Blue R-250, 25 g | GR135-25 |

Hoefer, Inc.

84 October Hill Road
Holliston, MA 01746

Toll Free: 1-800-227-4750

Phone: 1-508-893-8999

Fax: 1-508-893-0176

E-mail: support@hoeferinc.com

Web: www.hoeferinc.com

Chroma, Hoefer, and MiniVE are
trademarks of Hoefer, Inc.

Contrad 70 is a trademark of
Decon Laboratories.

Tween is a trademark of ICI
Americas Inc.

Coomassie is a trademark of ICI plc.

© 2012 Hoefer, Inc. —

All rights reserved.

Printed in the USA.